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** Easy Medium

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(0250S)

Medium packaged into stick pouch
For detection of coliform bacteria

** Easy Medium

BGLB Broth (Granule)

Product is divided powder media weighed and packaged into aluminum stick pouch for 400 mL. It can be used without weighing and dividing.

Formula (in 40.0g for 1 liter)

Calf Bile Powder	20.0g
Lactose	10.0g
Peptone	10.0g
Brilliant Green	0.0133g
	pH 7.0 ± 0.2

Directions

Put one pouch (16.0 g) into 400 mL of distilled water, mix well and heat to dissolve the medium. Distribute about 10 mL into individual medium size of test tube containing a fermentation tube or a Durham's tube. Sterilize by autoclaving at 121°C for 15 minutes, and cool down rapidly for use. Normally inoculate one (1) mL of test sample, and incubate it at 35 ± 1°C for 48 ± 3 hours.

When inoculate the test sample of more than 10 mL, use double concentration medium. (To prepare double concentration medium, put one pouch into 200 ml of distilled water.)

Determination and Counting

In case of examination of water, ice and snow, incubate in Lactose Broth first, and when the production of gas is observed, transplant the culture to BGLB Broth. When gas was evolved after incubation, coliform test is positive.

For the test of food and drinks, add the specimen directly to the medium and observe the production of gas.

Purpose

Product is used for detection of coliform bacteria in refreshing drink, milk and dairy products, meat and fish cake.

Warning and Direction for Use

1. General precautions

- 1) Read and follow precisely the warning and direction for use described on the package insert and/or label.
- 2) Do not use the product after its expiry date. Quality of the product is not warranted after its shelf life.

- 3) Do not use the product with damaged aluminum pouch or that contains any foreign materials.
- 4) After cut aluminum pouch, recommend using it at once. Avoid reuse after storage of prepared media.

2. Precautions for danger

- 1) When if medium touched eyes or mouth, immediately wash with plenty of water, and consult a physician.
- 2) Manipulations with microorganisms involve always certain risks of laboratory-acquired infections. Manipulations should be practiced under the supervision of key specialist with biohazard protection measures.

3. Precautions for disposal of waste

Any medium, materials and apparatus must be sterilized by autoclaving or boiling water after use, and then dispose as medical or industrial waste according to the Law on Waste Disposal and Cleaning. Also follow to local laws and regulations related to dispose such material after differentiation.

Storage

Keep dry at room temperature.

Shelf Life

Three (3) years after manufacturing. Actual expiry date is printed on label of box and aluminum pouch.

Package

Easy Medium BGLB Broth

For 400 mL (16.0g) x 60 pouches Code 56207

**/* Further information

Customer Support Section,

Shimadzu Diagnostics Corporation

3-24-6, Ueno, Taito-ku, Tokyo 110-0005 Japan

TEL: +81-3-5846-5707

E-mail: contact@sd.c.shimadzu.co.jp

** Manufactured by

Shimadzu Diagnostics Corporation

3-24-6, Ueno, Taito-ku, Tokyo 110-0005 Japan

** Revised : April 2023

* Revised : January 2018
(0250S)