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Medium packaged into stick pouch
For isolation of *Staphylococcus aureus*

** Easy Medium

Mannitol Salt Agar (Granule)

* Product is divided granulated media weighed and packaged into aluminum stick pouch for 200 mL. It can be used without weighing and dividing.

* Formula (in 111.0g for 1 liter)

This medium can be adjusted and/or supplemented according to the performance criteria required:

Beef Extract	1.0g
Peptone	10.0g
Sodium Chloride	75.0g
D-Mannitol	10.0g
Phenol Red	0.025g
Agar	15.0g
pH 7.4 ± 0.2	

* Directions

Open an aluminum pouch, put the medium (22.2 g) into 170 mL of distilled water, mix well and sterilize by autoclaving at 121°C for 15 minutes. Maintain the medium at about 50°C. Add 20% egg yolk solution to it and mix well. Distribute about 30mL of the medium into Petri dishes. Inoculate test sample, and incubate it at 37 ± 1°C for 48 ± 2 hours.

[Preparation of Egg Yolk Solution (ratio 20%)]

Fresh egg is sterilized by alcohol, put the egg yolk alone to sterilized cylinder aseptically. Add 80 mL of sterilized sodium chloride solution to the 20 g of egg yolk, and mix well to have homogeneous solution.

Determination and Counting

When staphylococci that utilize mannitol grow on the medium, they form yellow colonies and turn the medium around the colonies yellow, and so it is easily distinguished from bacteria that do not utilize mannitol. *Staphylococcus aureus* grown on Mannitol Salt Agar with Egg Yolk generate yellow (sometimes white) circle colony with pearl turbid ring (halo) around the colony by egg yolk reaction.

Purpose

Product is used for isolation of *Staphylococcus aureus* in foods or from environment of food production and cooking (apparatus, devices or workers fingers).

Staphylococcus aureus can be distinguished by capability of utilizing Mannitol and egg yolk reaction.

Warning and Direction for Use

1. General precautions

- 1) Read and follow precisely the warning and direction for use described on the package insert and/or label.
- 2) Do not use the product after its expiry date. Quality of the product is not warranted after its shelf life.
- 3) Do not use the product with damaged aluminum pouch or that contains any foreign materials.
- 4) After cut aluminum pouch, recommend using it at once. Avoid reuse after storage of prepared media.

2. Precautions for danger

- 1) When if medium touched eyes or mouth, immediately wash with plenty of water, and consult a physician.
- 2) Manipulations with microorganisms involve always certain risks of laboratory-acquired infections. Manipulations should be practiced under the supervision of key specialist with biohazard protection measures.

3. Precautions for disposal of waste

Any medium, materials and apparatus must be sterilized by autoclaving or boiling water after use, and then dispose as medical or industrial waste according to the Law on Waste Disposal and Cleaning. Also follow to local laws and regulations related to dispose such material after differentiation.

* Storage

Keep dry at room temperature (1~30°C).

Shelf Life

Three (3) years after manufacturing. Actual expiry date is printed on label of box and aluminum pouch.

** Package

Easy Medium Mannitol Salt Agar
for 200 mL (22.2g) x 46 pouches

Code 56213

** Further information

Customer Support Section,
Shimadzu Diagnostics Corporation
3-24-6, Ueno, Taito-ku, Tokyo 110-0005 Japan
TEL: +81-3-5846-5707
E-mail: contact@sd.c.shimadzu.co.jp

** Manufactured by

Shimadzu Diagnostics Corporation

3-24-6, Ueno, Taito-ku, Tokyo 110-0005 Japan

** Revised : April 2023

* Revised : August 2020
(0H01S)